DOC272.98.80487

**97-0073-02, 97-0073-03,  
98-0002-01, 90-0012-01,  
99-0064-03, 90-0086-03**

09/2014, Edition 1

**Safety Information User Instructions  
Sicherheitsinformationen/Anweisungen für Benutzer  
Informazioni di sicurezza - Istruzioni per l'utente  
Informations relatives à la sécurité et instructions pour l'utilisateur  
Instrucciones de uso sobre información de seguridad  
Instruções do utilizador sobre informações de segurança  
Gebruikershandleiding veiligheidsinformatie  
Sikkerhedsoplysninger brugerinstruktioner  
Informacje dotyczące bezpieczeństwa – instrukcje dla użytkowników  
Säkerhetsinformation användarinstruktioner  
Turvaohjeet käyttäjälle  
Biztonsági előírások felhasználóknak  
Instrucțiuni de siguranță pentru utilizator  
Saugos informacija ir instrukcijos naudotojams  
Указания по безопасности — инструкции для пользователя  
Güvenlik Bilgileri Kullanıcı Talimatları  
Bezpečnostné informácie – návod na použitie  
Varnostni napotki za uporabnike  
Οδηγίες χρήστη και πληροφορίες για την ασφάλεια  
Informacije o bezbednosti za korisnike**

English .....	3
Deutsch .....	5
Italiano .....	7
Français .....	9
Español .....	12
Português .....	14
Nederlands .....	17
Dansk .....	19
Polski .....	21
Svenska .....	23
Suomi .....	25
Magyar .....	27
Română .....	29
lietuvių kalba .....	31
Русский .....	33
Türkçe .....	36
Slovenský jazyk .....	38
Slovenski .....	40
Ελληνικά .....	42
Српски .....	45

# USER INFORMATION FOR PEN-RAY LAMPS

**IMPORTANT NOTICE:** When using a Pen-Ray® light source, the following instructions are recommended to insure proper lamp operation.

## WARNING

Exposure to UV light can cause eye and skin damage. Wear appropriate protective eyewear equipment when the lamp is on.

1. Use the recommended power voltage and current.
2. Consult UVP for feasibility of any unusual applications.
3. Good air circulation around the lamp is recommended.
4. Do not inhibit air circulation around handle.
5. Consult UVP regarding any extended lead cord lengths.
6. **CAUTION** : Pen-Ray Lamps produce UV radiation which is harmful to unprotected eyes and skin. Protective glasses and clothing are recommended.
7. Quartz material is fragile - HANDLE WITH CARE
8. If this product contains Mercury (see box label), handle with care and dispose of properly. MSDS available upon request.

**WARRANTY:** Pen-Ray Lamps are warranted for 90 days from date of purchase. Contact UVP for information on Pen-Ray power supplies, UV intensity meters and UV blocking eyewear.

## PEN-RAY POWER SUPPLIES

PS-1, PS-2, PS-3, PS-4, PS-9 and PS-11

## WARNING

Please read operating instructions before turning unit on.

## INTRODUCTION

The Pen-Ray Power Supplies are made specifically for the Pen-Ray Lamps. These power supplies provide proper voltage and current for starting and operating the lamps. Mating recommendations are available to provide optimum energy output from and maximum life span of the Pen-Ray lamps.

## DESIGN

Each power supply is designed to be maintenance free. There is an internal circuit protection fuse which is designed to prevent mains circuit shut down in the unlikely case of a catastrophic internal failure. The connector must be used with the mating connector supplied with the Pen-Ray lamps. The connector is certified to **UL** and **CE** compliance for use with this power supply. The power supply carries the **CE**, **ETL** and **ETLc** marks.

## OPERATING INSTRUCTIONS

The operation of the PS series power supply is as follows:

1. Make sure the switch is in the off position.
2. Connect the Pen-Ray lamp to the connector making sure the connector is completely inserted into its mating part.
3. Plug the unit into proper mains power supply.
4. Turn on the unit and the lamp should immediately light up.
5. Always unplug the unit when disconnecting or connecting the lamp.

## MAINTENANCE

Keep the connector clean and free from debris, oil, fluids, etc. If the cord should become frayed or cracked, it should be replaced immediately. Call UVP for replacement service.

If connector arcing should occur due to dirt in the connector or attempting to operate a Pen-Ray lamp which operates in excess of 300 volts, do not continue operation. The unit should be returned to the factory for repair. (Cleaning the carbonized area between the connector pins may restore function to the power supply).

## WARNINGS

- Never handle the connector when the unit is in position as a hazardous shock condition will exist.
- Do not attempt to operate lamps with operating voltages greater than 300 volts continuously. Starting voltages do not apply.
- If the equipment is used in a manner not specified by UVP, LLC, the protection provided by the equipment may be impaired.

### Unit may become HOT during continuous use.

If internal circuit protection fuse blows, the unit should be returned to the factory for service. The internal protection fuse is designed to trip if the primary coil short circuits for any reason, protecting the user and the circuit it is plugged into. The unit may be cleaned with any mild cleaning agent. Although the powder coated finish is extremely durable, avoid acids and caustic substances. Do not use solvent cleaners on plastic end caps.

## ENVIRONMENTAL CONDITIONS

- For indoor use only
- Temperature: 5 to 40°C
- Maximum or Relative Humidity: 80% for temperatures up to 31°C decreasing linearly to 50% relative humidity at 40°C
- Mains supply voltage fluctuations not to exceed +10% of the nominal voltage
- Installation is category II
- Pollution degree is 2

## ULTRAVIOLET BLOCKING EYEWEAR

**Important:** UV Eyewear is intended to protect the eyes from harmful ultraviolet radiation emitted from UV lamps. The Eyewear is designed for eye protection from lamps emitting wavelengths of 254nm to 365nm (nanometers). These UV Blocking Eyewear also reduce longwave ultraviolet "blue haze" for optimum viewing.

**Caution:** Although the Goggles and Spectacles are intended to protect the eyes and the Faceshield to protect the face and neck, care must be taken to protect other parts of the body from skin burn associated with shortwave UV wavelengths. Avoid shining any ultraviolet lamp on reflective surfaces. Wear protective clothing and skin cream.

Eyewear meet performance standards of ANSI Z87.1-1989. Spectacle and Goggle lenses are impact resistant, but not unbreakable. Clean all eyewear surfaces with a mild soap solution and inspect frequently. Do not use solvent on lenses. Carefully rinse off all soap. Air dry or pat dry with a clean, soft cloth or tissue. Pitted or scratched eyewear reduce protection. If scratched or broken, replace immediately.

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P.O. Box 389, Loveland, CO 80539-0389 U.S.A.
(800) 227-4224 (U.S.A. only)